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**Sample Report**
Result Manager

<div style="border: 1px solid black; padding: 2px;"> <b>Contact Information</b>                  Emergency Room                  Hospital Name                  123 Main Street                  Anywhere DC 123456                  USA             </div>	<div style="border: 1px solid black; padding: 2px;"> <b>Dr Generic Contact Information</b>                  Phone: 555-555-5555             </div>	[ LOGO ]																																																						
<b>Patient Information</b> Patient ID: 200 Name: Smith, John R      Age (y): 86      Height (cm): 62 Sex: M      Date of Birth: Jun 8, 1928      Weight (kg): 100																																																								
<b>Sample Information</b> Sample ID 100 Volume 100.00 mL Sample Type S																																																								
<table border="1" style="width: 100%; border-collapse: collapse;"> <tr> <td style="width: 50%;">Release Date/Time</td> <td>Sep 9, 2014 4:37:25 PM</td> </tr> <tr> <td>Collected Date/Time</td> <td>Sep 9, 2014</td> </tr> </table>			Release Date/Time	Sep 9, 2014 4:37:25 PM	Collected Date/Time	Sep 9, 2014																																																		
Release Date/Time	Sep 9, 2014 4:37:25 PM																																																							
Collected Date/Time	Sep 9, 2014																																																							
<table border="1" style="width: 100%; border-collapse: collapse;"> <thead> <tr> <th>Test</th> <th>Aspect</th> <th>Result</th> <th>Unit</th> <th>Prio.</th> <th>Reference Range</th> </tr> </thead> <tbody> <tr> <td>Acetaminophen (ACTM)</td> <td></td> <td>110</td> <td>ug/mL</td> <td>S</td> <td>[ 10.0 - 30.0 ]</td> </tr> <tr> <td colspan="6"> <i>Result Comment:</i> Instrument Result Comment                      EasyLink Result Comment                 </td> </tr> <tr> <td>Thyroid Stimulating Hormone (TSH)</td> <td>DOSE</td> <td>50</td> <td>ng/mL</td> <td>R</td> <td>[ - ]</td> </tr> <tr> <td colspan="6"> <i>Result Comment:</i> Centaur User Comment                 </td> </tr> <tr> <td>Thyroid Stimulating Hormone (TSH)</td> <td>COFF</td> <td>1.0</td> <td>ng/mL</td> <td>R</td> <td>[ - ]</td> </tr> <tr> <td colspan="6"> <i>Result Comment:</i> Centaur User Comment                 </td> </tr> <tr> <td>Thyroid Stimulating Hormone (TSH)</td> <td>RLU</td> <td>4482</td> <td></td> <td>R</td> <td>[ - ]</td> </tr> <tr> <td colspan="6"> <i>Result Comment:</i> Centaur User Comment                 </td> </tr> </tbody> </table>			Test	Aspect	Result	Unit	Prio.	Reference Range	Acetaminophen (ACTM)		110	ug/mL	S	[ 10.0 - 30.0 ]	<i>Result Comment:</i> Instrument Result Comment EasyLink Result Comment						Thyroid Stimulating Hormone (TSH)	DOSE	50	ng/mL	R	[ - ]	<i>Result Comment:</i> Centaur User Comment						Thyroid Stimulating Hormone (TSH)	COFF	1.0	ng/mL	R	[ - ]	<i>Result Comment:</i> Centaur User Comment						Thyroid Stimulating Hormone (TSH)	RLU	4482		R	[ - ]	<i>Result Comment:</i> Centaur User Comment					
Test	Aspect	Result	Unit	Prio.	Reference Range																																																			
Acetaminophen (ACTM)		110	ug/mL	S	[ 10.0 - 30.0 ]																																																			
<i>Result Comment:</i> Instrument Result Comment EasyLink Result Comment																																																								
Thyroid Stimulating Hormone (TSH)	DOSE	50	ng/mL	R	[ - ]																																																			
<i>Result Comment:</i> Centaur User Comment																																																								
Thyroid Stimulating Hormone (TSH)	COFF	1.0	ng/mL	R	[ - ]																																																			
<i>Result Comment:</i> Centaur User Comment																																																								
Thyroid Stimulating Hormone (TSH)	RLU	4482		R	[ - ]																																																			
<i>Result Comment:</i> Centaur User Comment																																																								
<b>Sample Comments:</b> Instrument Sample Comment																																																								
Sep 9, 2014 _____ Signature																																																								

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## Sezione 2: User-Initiated Sample and Result Action Buttons

**Table 3. User-Initiated Sample and Result Action Buttons: Location and Description**

Action Button	EasyLink User Interface Page	EasyLink User Interface Page Section	Button Name
Release Patient	Patient	All	Release
QC Result	QC Today's Result	All	Release
Release Sample	Sample Review	Sample Details	Release
Release Result	Sample Review	Results	Release
Release Result	Result Details	Result Value	Release
Release Result	Sample Details	Results	Release
Release All Results	Sample Details	Results	Release All
Delete Result	Sample Review	All	Delete
Reset Result	Sample Review	All	Reset
Reorder Result	Sample Review	All	Reorder
Ignore Result	Sample Review	All	Ignore

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Figura 2. EasyLink User Interface Sample and Result Action Button Details

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1. Sample Details section	7. Ignore Result button
2. Release Sample button	8. Results section
3. Delete Result button	9. Results scrollbar*
4. Release Result button	10. Results Measurement Time column
5. Reorder Result button	11. Results Status column
6. Reset Result button	12. Results Checkbox column

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### **Limited Use of User-initiated Sample and Result Action Buttons**

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**Siemens Healthcare Diagnostics S.r.l.**